## **Product Datasheet**

# Lightning-Link (R) Rapid Type A Biotin Antibody Labeling Kit 370-0030

Unit Size: 3 x 10ug Reaction

Store at -20C.

www.novusbio.com



technical@novusbio.com

**Publications: 3** 

Protocols, Publications, Related Products, Reviews, Research Tools and Images at: www.novusbio.com/370-0030

Updated 10/23/2024 v.20.1

Earn rewards for product reviews and publications.

Submit a publication at www.novusbio.com/publications Submit a review at www.novusbio.com/reviews/destination/370-0030



#### 370-0030

Lightning-Link (R) Rapid Type A Biotin Antibody Labeling Kit

Product Information	
Unit Size	3 x 10ug Reaction
Concentration	Concentration is not relevant for this product. Please see the protocols for proper use of this product.
Storage	Store at -20C.
Conjugate	Biotin
Buffer	

#### **Product Description**

#### **Description**

Lightning-Link Rapid Antibody Labeling kits enable direct labeling of antibodies, proteins, peptides or other biomolecules (See protocol for further information) and allow conjugations to be set up in seconds, and used within minutes (less than 20 minutes).

Our Lightning-Link Rapid technology can be used to label small quantities of antibody with 100% recovery. The researcher simply pipettes the antibody or other biomolecule into a vial of lyophilized mixture containing the label of interest and incubates the reaction for **only 15 minutes** (for more details please watch the video below). As there are no desalting or dialysis steps that commonly interrupt traditional protein labeling procedures.

**Features Benefits** Quick and easy to use Save time, no special knowledge required No separation steps 100% recovery - no antibody/protein loss Can be used in a wide range of applications Flexible Freeze dried Ships at ambient temperature, long shelf-life Fully scalable (10 ug to 1 g or more) Easy transfer from R&D to manufacturing Stringently QC tested Consistent high quality, excellent batch-to-batch reproducibility Large number of labels available Experimental flexibility Reliable: nearly 300 references Successfully used in many fields of research

The attachment of biotin to biomolecules is an important laboratory technique. Biotin binds to the tetrameric avidin proteins, including streptavidin and neutravidin, with exceptionally high affinity, and this interaction is exploited in various applications such as western blotting, immunohisthochemistry and ELISA. Our biotinylation kits are optimised to create high quality conjugates and minimise assay background.

#### Lightning-Link Biotin Has been optimized for two separate applications:

- Type A Intended for assays in which a streptavidin-labeled detection reagent will be used.
- Type B Optimised for assays in which the biotinylated protein is captured by streptavidin immobilized on a surface (ie plates, nitrocellulose, magnetic beads etc).

Learn more about Lightning-Link™ Conjugation Kits by reading FAQs

For more information please check out these useful links!

<u>Antibody Labeling Guide</u>

<u>Antibody Conjugation Illustrated Assay</u>




Kit Components	1 or 3 glass vial(s) of Lightning-Link Rapid mix, 1 vial of LL-Rapid Modifier reagent, 1 vial of LL-Rapid Quencher reagent
Notes	This product is manufactured by Abcam and distributed by Novus Biologicals.  This product is for research use only and is not approved for use in humans or in clinical diagnosis. This product is guaranteed for 1 year from date of receipt and this statement overrides any mentioned guarantee period on the limitations section of this products datasheet. Please contact technical@novusbio.com with questions.
<b>Product Application Details</b>	
Application Notes	By circumventing the desalting or dialysis steps that commonly interrupt traditional antibody conjugation procedures, LightningLink technology can be used to label both small (e.g. 10 ug) and large quantities of primary antibodies

suitable for labeling up to 20 ug of antibody.

with ease. Batch-to-batch variation upon scale up is minimal as the process is so simple, and recoveries are always 100%. This kit is supplied with 3 vials, each

#### **Images**

Lightning-Link Rapid Type A Biotin Antibody Labeling Kit [370-0030] - CRP labeled with Lightning-Link Rapid Biotin A at 10ug, 10ug, 1mg, 10mg and 100mg single reaction vials tested with Streptavidin-Rabbit IgG in a sandwich ELISA. Detection: ABTS.

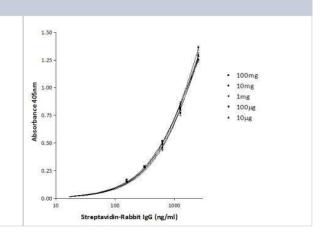

#### **Publications**

Hughes AC, Kirkland M, Du W et al. Development of Thermally Stable Nanobodies for Detection and Neutralization of Staphylococcal Enterotoxin B Toxins 2023-06-16 [PMID: 37368700]

Noreng S, Ota N, Sun Y et al. Structure of the core human NADPH oxidase NOX2 Nature communications 2022-10-14 [PMID: 36241643]

Laoharawee K, DeKelver RC, Podetz-Pedersen KM et al. Dose-Dependent Prevention of Metabolic and Neurologic Disease in Murine MPS II by ZFN-Mediated In Vivo Genome Editing. Mol. Ther. 2018-03-10 [PMID: 29580682] (WB)





### Novus Biologicals USA

10730 E. Briarwood Avenue Centennial, CO 80112 USA

Phone: 303.730.1950 Toll Free: 1.888.506.6887

Fax: 303.730.1966

nb-customerservice@bio-techne.com

#### **Bio-Techne Canada**

21 Canmotor Ave Toronto, ON M8Z 4E6 Canada

Phone: 905.827.6400 Toll Free: 855.668.8722 Fax: 905.827.6402

canada.inquires@bio-techne.com

#### **Bio-Techne Ltd**

19 Barton Lane Abingdon Science Park Abingdon, OX14 3NB, United Kingdom Phone: (44) (0) 1235 529449

Free Phone: 0800 37 34 15 Fax: (44) (0) 1235 533420 info.EMEA@bio-techne.com

#### **General Contact Information**

www.novusbio.com Technical Support: nb-technical@biotechne.com

Orders: nb-customerservice@bio-techne.com

General: novus@novusbio.com

#### Limitations

This product is for research use only and is not approved for use in humans or in clinical diagnosis. Kits are guaranteed for 6 months from date of receipt.

For more information on our 100% guarantee, please visit www.novusbio.com/guarantee

Earn gift cards/discounts by submitting a review: www.novusbio.com/reviews/submit/370-0030

Earn gift cards/discounts by submitting a publication using this product: www.novusbio.com/publications

